# DigiMicro Mobile

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<b>Security instructions</b> .....	29
<b>Appliance description</b>	
System requirements .....	30
Delivery content .....	30
Appliance description.....	31
Getting started .....	32
Charging the battery .....	33
Insert the SD card .....	34
Photo mode.....	35
Turn on/off .....	35
How to capture .....	36
Record video .....	37
Playback.....	38
Picture settings .....	40
Basic settings .....	42
Software installation .....	44
Picture editing .....	47
Additional information .....	49
Magnification .....	49
Mounting the stand .....	50
Service informationen .....	50
Specifications .....	51
Declaration of conformity .....	51
Disposal of electronic equipment .....	51
Warranty regulations .....	103

## The company placing the product on the market

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**Please read this information carefully before you use the appliance.**

### **General instructions**

- Don't place the unit on unstable location, if it falls down, people may be injured.
- The unit is not a toy, please beware of small children.
- If you connect the unit to other electronic devices, please study also the security instruction of this device.
- If you have any question regarding the unit, how it works, the safety or the correct connection please contact our technical support or ask any specialist.
- The seller will not be liable for any damages caused by misuse of the product and any claim from third parties.

### **Environmental conditions**

- Don't drop or shake the unit, it might be damaged.
- Avoid too high pressure to the LC-display, it might be damaged.
- Don't leave the unit where the temperature is higher than 60°C, e.g. in cars with closed windows in the summertime, near heaters or any other type of heat source.
- Don't use the unit in rooms with high humidity e.g. bathrooms.
- Don't use the unit at dusty places.

### **Technical problems**

- In case any foreign material or liquid gets into the product, please immediately remove the power cord. Before you use it again, ask your authorised dealer to check the product carefully.
- Please do not hesitate to contact our service center.

### **Maintenance**

- We strongly suggest using a clean, dry, non-alcohol cotton swab to clean the glass surface periodically to produce a clear image during operation.

### **Batteries**

- If a battery has leaked, avoid any contact with your skin, eyes or mucous membranes. If necessary, rinse the affected parts with water and consult a doctor immediately.

### **DANGER to children**

- Children are not aware of the risks involved due to improper handling of electronic appliances. Therefore, keep children away from the product.
- Batteries can be highly dangerous if swallowed. For this reason, keep batteries and remote control out of the reach of small children. If a battery has been swallowed, seek immediate medical advice.
- Keep children away from the packing material. There is a risk of suffocation!

## Delivery content

### System Requirements

- OS: MS Windows XP SP2/Vista/ Windows7 and MAC OS ab 10.6 (without measuring software)
- min. processor Pentium 1800 or compatible AMD processor
- 100 MB free hard disk space
- min. 32MB RAM, 512 MB RAM recommended
- 24 Bit colour Video adaptor
- free USB 2.0 port
- CD-ROM drive

### Delivery content

DigiMicro Mobile

stand

charger

AV cable

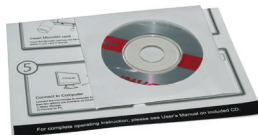
USB cable

Carry bag

user manual


CD and QG


Cleaning fabric

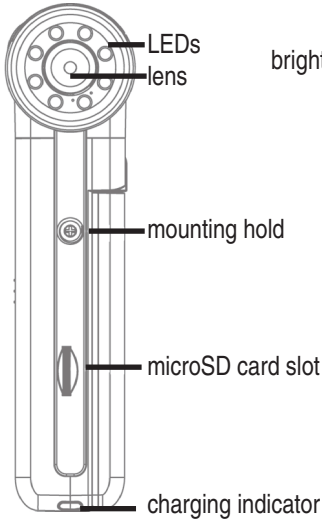
*View of the device*


*front*

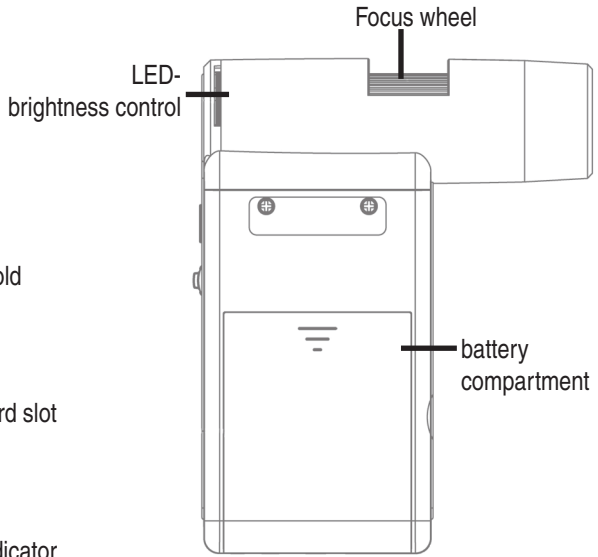*right side*

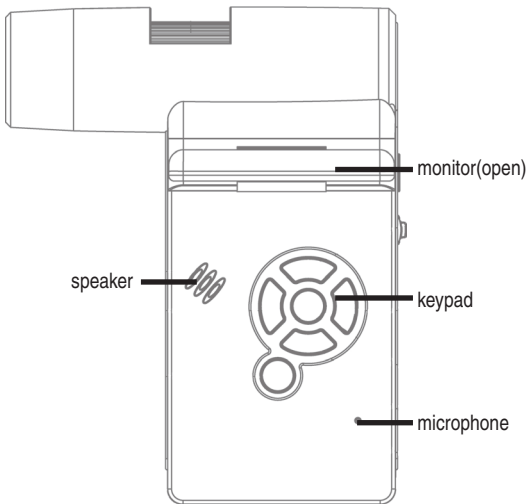
*left side*

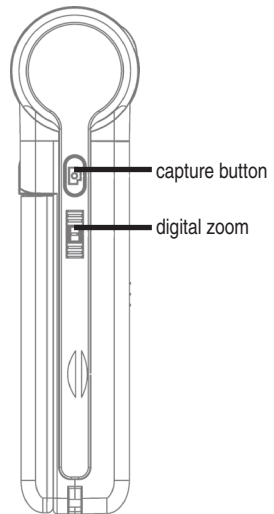*rear side*


## View of the device
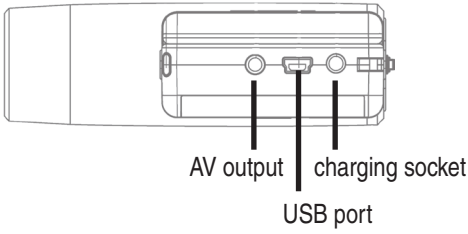
### View of the device


**bottom**
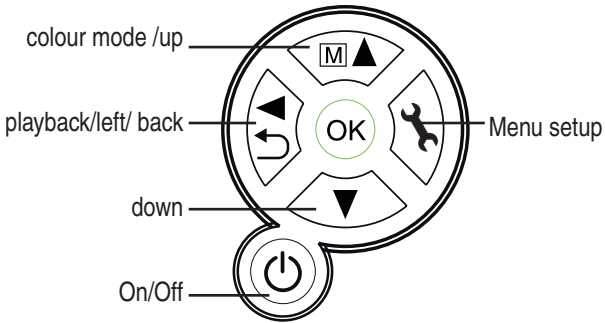
### Keypad and navigation

### Getting started
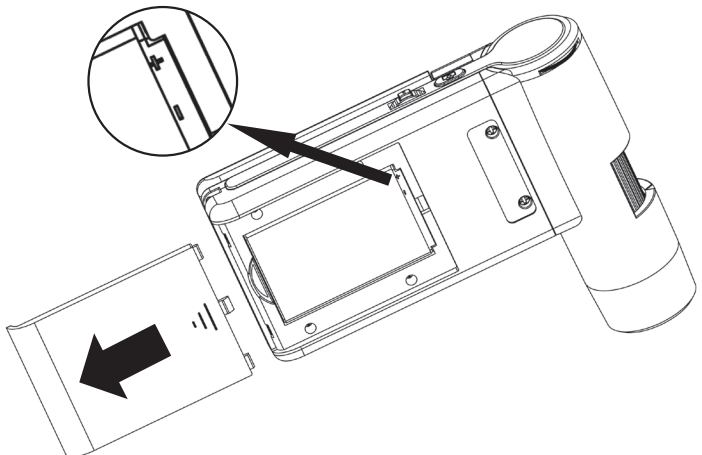
#### Install the battery


Open the battery door by sliding it outward. Insert the lithium battery to the battery compartment according to the polarities ⊕ ⊖ shown in the compartment.

### Charging the battery

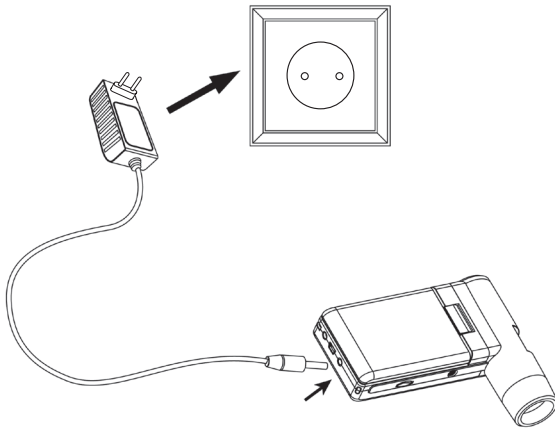Before using the microscope, you must charge its battery.

- Plug the small end of the power adaptor into the DC-in port located on the bottom side of the unit. Then plug the other end of the power adaptor into a standard AC outlet.
- The battery indicator LED at the end corner of the unit will be in red to show that the battery is charging.
- When the battery is finished charging, the battery indicator LED will be in off. Unplug the unit.
- You can now use your microscope for up to four hours of continuous use. When the battery is low, there will be low-power indicator on the right down corner of the panel, front LED light will become weak and image quality becomes poor. Then you have to connect the power adaptor and plug it in to recharge the battery.

### Battery charging indicator

The display shows the charging level on the right side. Additional indicators are the LED brightness and the image quality. If the battery is empty, the brightness goes down and also the image quality.

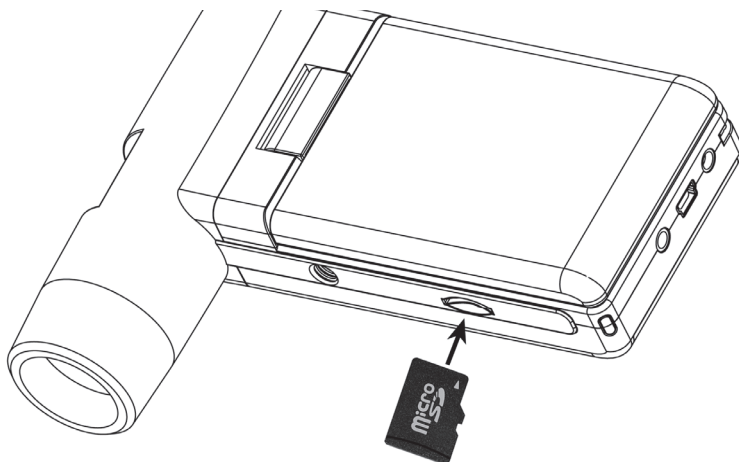
### Insert a MicroSD Card and the SD card indicator

You must insert a MicroSD card to record and store your pictures and videos.

- Turn the device off.
- Insert the MicroSD card into the MicroSD card slot; gently press in until card is latched.
- The display shows a SD card icon.


Please format the card before the first usage. **(Format SD)**.

**Do not force the card into the slot; forcing may damage microscope and the MicroSD memory card. Align MicroSD card position (as marked) and try again. The MicroSD memory card must be formatted before capture.**

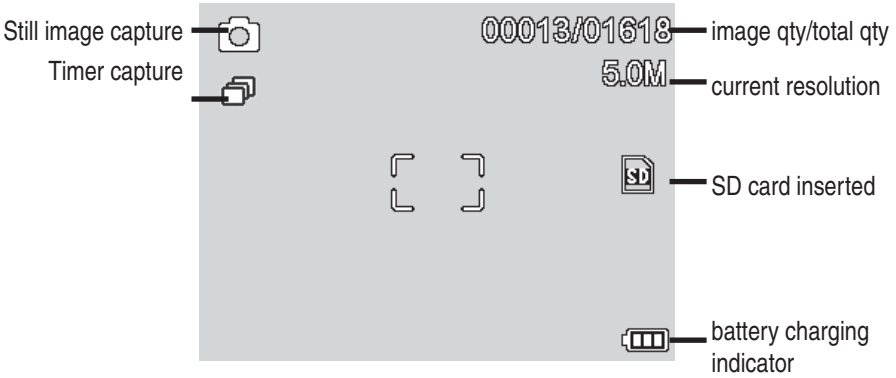
**Display - photo mode**

The icon displays can be switched on/off by pressing Zoom Wheel downward vertically.

**Turn on**

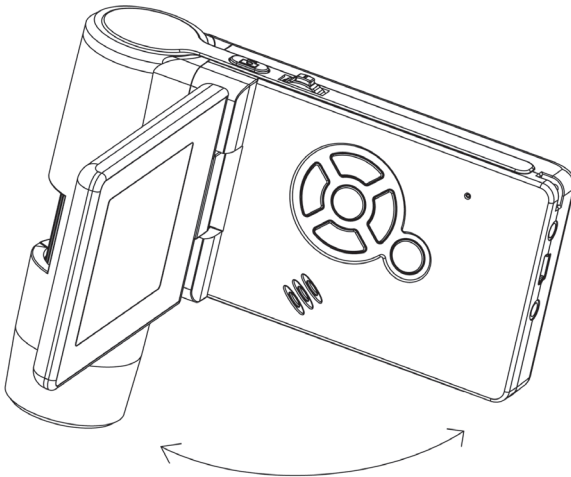- Unfold the display panel, and then the microscope will power on directly.
- If need to power on when panel unfolded, press **POWER ON/OFF** button.


**Turn off**

- Fold the display panel and then the microscope will turn off automatically
- Press ON/OFF button
- By Power Save.


## Using the microscope

### How to capture

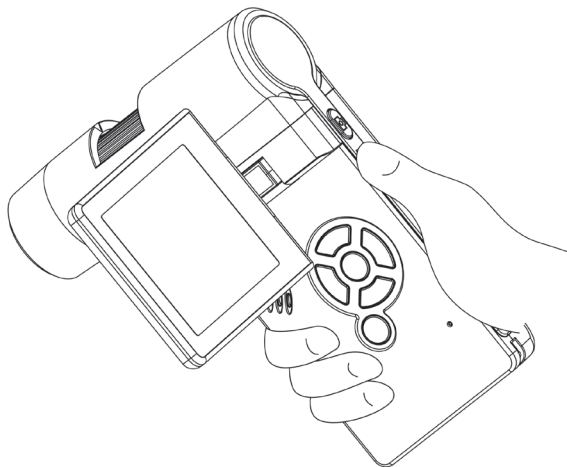
- Turn on the microscope by simply unfolding its display panel to a proper position, normally at an angle of 90 degree with the viewing end as follows:

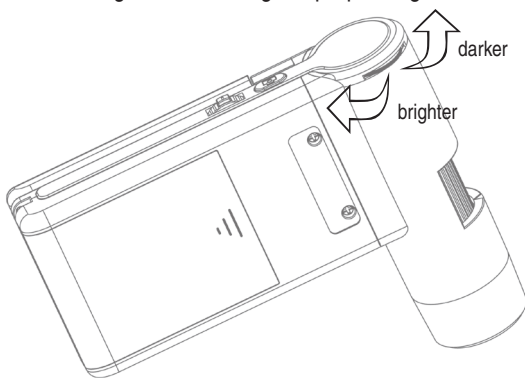- Hold the microscope main body by right hand; put the viewing end closely to target object.
- Adjust the LED brightness dial to get a proper brightness for the LEDs at the viewing end.

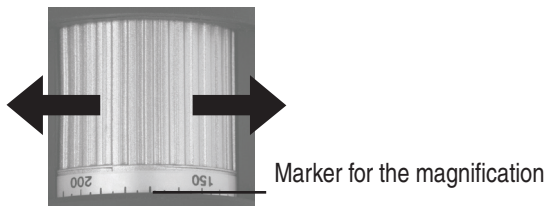
- Adjust focus wheel to get clear image by left hand.

- When a clear image got, hold the microscope steadily and press Capture button one time.

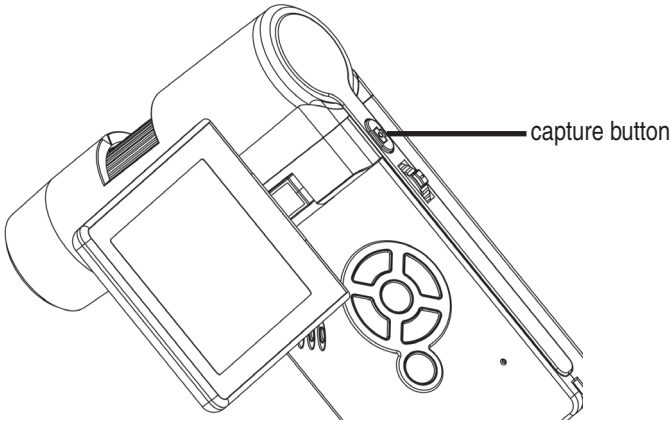- Press the capture button

**i** When a clear image got, hold the microscope steadily and press Capture button one time. Please see timer on page 40 also.  
Remember the magnification on the scale for later measurements.

**Record Video**

Press ◀ / ▶ button to enter video mode as follows. There is video icon  on the left corner.

- Press **Capture button** to start recording video.
- Press again to stop.

- The video has 2 resolution: VGA (640x480 pixel) and QVGA (320x240 pixel).
- To change the resolution press  button and select VGA or QVGA. To leave press the **button** again.

### Playback

To enter the playback mode press the -**button** until the display shows the playback icon . In case the SD card contains video files only, the display shows the video icon .

- To view images and videos press the **buttons**.
  - The microscope is able to recognize the file type, picture or video, and shows an icon accordingly.

- The current selected file is a picture.
- Select the previous and next picture with the **buttons**.

- The current selected file is an video.

### Navigation:

**OK**:▶ Start the playback with the **OK button**  
For pause press **OK button** again.

**fast forward and rewind**: For fast forward and rewind press the **buttons**. To switch back to normal playback press the **OK button**. If the last choice was „rewind“ it plays in a normal speed but backward. To switch to normal playback (forward) press once the forward button and **OK**.

**OK**:▶ Finish the fast forward and rewind.

### Playback menu

During the playback you will get access to a submenu with the following functions.

Delete, protect and slide show. To access the playback menu press the **button**.

### Navigation:

**buttons** : Select a menu item

**button** : Exit the menu.

**OK button** : Confirm the selection.

### Delete:

In the following submenu select

**This image:** delete the image shown in the display.

**All:** Delete all

To delete confirm the prompt with **OK(Execute)**.

**Protect for image and video files.**

Protect/ Unprotect pictures and videos from delete.

**Lock one:** Protect the picture in the display.

**Unlockm One:** Remove protection of the picture in the display.

**Lock All:** Protect all pictures

**Unlock all:** Remove protection from all pictures.

**If a picture is protected the display shows a key icon.**

**Slide show:**

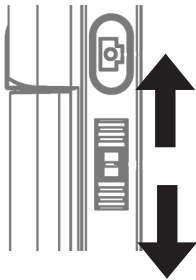
This menu item starts a slide show. A submenu shows the available intervals. To quit press the **OK button**.

**Colour mode** (only available in photo mode)

At preview status, press **[M]▲ button** to switch among Colour, Positive and Negative.

**Zoom**

The zoom wheel is located on the rear. The zoom range is 1.0 to 4.0 times. The zoom factor is shown in the display on the left side.

Zoom in, push Zoom wheel upward

Zoom out, push Zoom Wheel downward

**In playback mode the display shows thumbnails of the pictures when the zoom wheel is pushed downwardly.**

**Video records:** Push the zoom wheel downward if the playback is abnormal terminated (Pause) or you want to finish the playback. In this case you will return to the overview of the pictures. If the playback is finished you can select the previous and next file with the **▼/▲** buttons.

## Using the microscope

### Main menu - picture settings

Press the  **button** to enter the main menu.

Navigation:

 **buttons** : Select a menu item.

 **button** : In menu item back

 **button** : Exit the menu.

 **button** : Confirm the selection.

#### Resolution

The available resolution is VGA, 1.3, 3.0, 5.0, 9.0 and 12 megapixel.

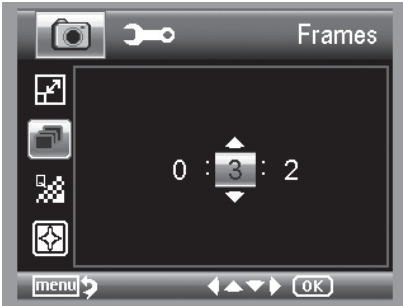
#### Timer

From the main menu, choose Timer and then enter Timer setup:

- **Frames**

Input the quantity of frames you want to take by ▼/▲ **buttons**

The above setup means 32 frames.

- **Intervall**

Input the interval time between frame capture by ▼/▲ **buttons**.

### Quality

The three selections will result in different image compress ratio.

### Sharpness

Choose Sharpness and then select among Super Fine/Fine/Normal.

## Using the microscope

### Main menu - basic settings

Press the  **button** to enter the main menu. drücken Sie die  **button**. Press the  /  **button** to enter the basic settings.

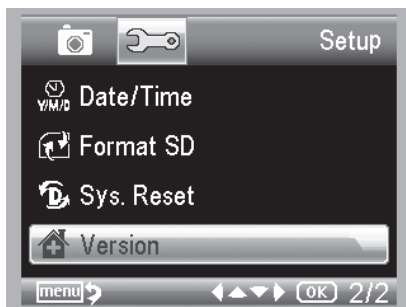
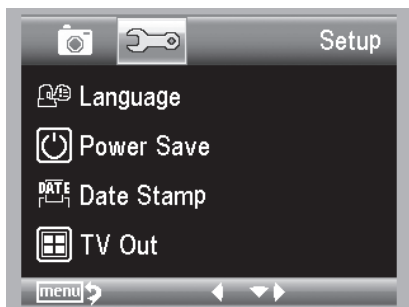
 /  **buttons** : Select a menu item.

 **button** : In menu item back

 **button** : Exit the menu.

 **button** : Confirm the selection.

English

### Language

Choose Language

### Power save

Choose Power Save and then select among the following a target time for the microscope to power off without key operation.

**Date Stamp**

Date/Time: both date and time will be marked on images taken.

Date: on date will be marked on images taken

Off: None of date and time info will be marked on images taken.

**TV Out**

Choose TV Out and then select, between NTSC and PAL, the correct mode compatible with your TV.

**Date/time**

Choose Date/Time to setup date and time.

Setup date and time by **▼/▲, ◀/▶ ↻-buttons.**

Setup date mode among YY/MM/DD, DD/MM/YY, and MM/DD/YY with **◀/▶ ↻-button.**

**Format SD card**

Choose Format SD to format the SD card before the first usage. Confirm the prompt with **OK(Execute)** otherwise select **Cancel**.

## Software installation

### Default settings

Choose **Sys. Reset** to reset the system back to default..

### Version

Choose Version to view the current machine version.

### Software installation

 **Install the software first before the microscope is connected to the USB port.**

- Insert the driver CD into CD-ROM Drive and this will automatically display the following interface:

- Click **Install ProtatableCapture** software and then reboot the system when asked.
- You can browse the User's manual in PDF format and the CD contents by choosing the corresponding menus.

## Download to Computer

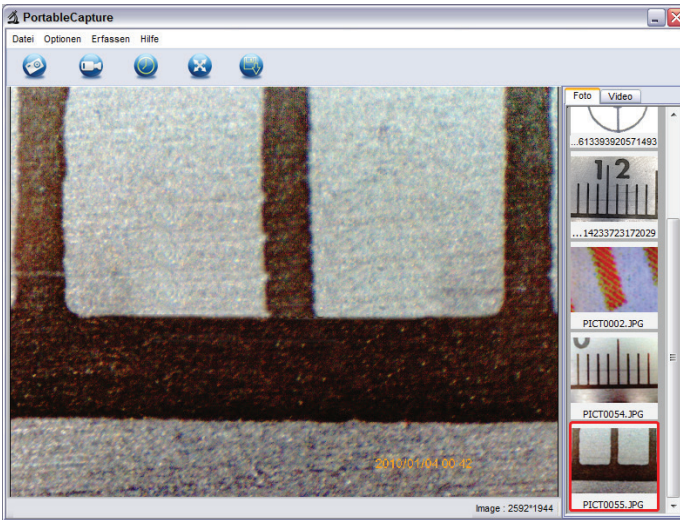
Connect the microscope to computer by USB cable. Choose Mass Storage and then you can view stored photos & videos.

Unfold the display panel. When working on computer, the button functions are no longer valid, but most of functions are available by PortableCapture software.

## Preview on Computer

- Connect the microscope to computer by USB cable.
- Choose PC Camera mode.
- Start PortableCapture software.

- When the Microscope is disconnected from your PC USB port, the screen will display “No Device detected, please connect your Microscope directly to your PC USB port.”

- If the mouse cursor is over a thumbnail on the right side the images is shown in the preview window. Click the right mouse button to enter a submenu where you can open, store copy an delete an image. During the storing procedure the program shows a save file dialog. The file can be saved in either jpeg or bmp. The jpeg size in which the photo is saved is the maximum allowed for jpeg. You can choose to decrease the size by adjusting quality. Click the Advanced button to view and select jpeg save options.

**In the PC camera mode the image size is limited to 320x240 and 640x480 pixel.**

- To capture a picture press F11 or clock the camera icon. The photos are stored in the **Photos Directory**.
- The images appears in the image list on the right side.
- To start record a video click on the video camera icon on the top of the window or select **Capture > Video**.
- To stop the record click once again on the video camera icon.
- The videos are stored in the **Videos Directory**.

### Main menu

#### File

- |                   |  |
|-------------------|--|
| Photos Directory: | Directory to store the photos.   |
| Videos Directory: | Directory to store the videos.   |
| Import:           | Import the pictures from the microscope to the computer for further use. |
| Exit:             | Finish the program.  |

#### Options

- |                    |   |
|--------------------|---|
| Preview size:      | In this menu you can set the preview size.  |
| Rotate             | Here you can rotate the image 90 degrees clockwise or   |
| Date/Time:         | The date and time is displayed in the image   |
| Language:          | Sets the OSD language.  |
| Toggle Fullscreen: | Switching between normal and screen display. The switch back to normal view is done with a double click on the image display.   |
| Preferences:       | Here you can make different settings for image display and different effects can be produced (only VGA resolution and smaller). The settings can be loaded and saved. The words on the screen are self-explanatory. |

#### Capture

- |        |  |
|--------|--|
| Photo: | Capture a single image                                   |
| Video: | Starts a video recording. Click again to stop recording. |

**As soon as the program is started and the microscope is not connected, an error message appears**

**“No Device detected, please connect your Microscope directly to your USB Port.”**

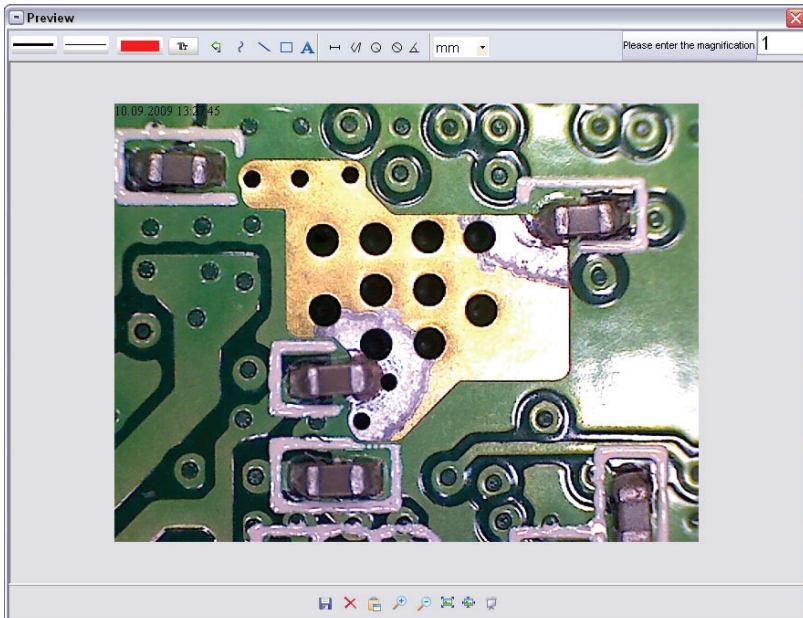
If your computer has no USB 2.0 interface, the image resolution decreases.

### Picture editing

Simultaneously with the main program also a software is installed, that allows you to perform measurements in the picture. To start the software, double click on the image in the picture gallery on the right side.

At the bottom of the window you see some icons, which you remember from the fax / Bildbe-trachter Windows. You can save the edited picture, delete, copy into the memory, increase or decrease the size.

Please enter the magnification

1

First set the magnification, which is read on the scale on the knob. On the cabinet there is a small indication. Simply read the value which points to the indication

## Line type

Set the desired line type

## Line width

Set the line width

## Line Color

## Character set for the dimension

## Cancel last edition

## Freehand drawing

## Draw a line (with no dimension)

## Draw a Rectangle

## Text entry mode

Distance measured with an indication of the distance between 2 points

## Multiple-path measurement

Radius Circle, measures the distance from the center of the circle to its radius.

## Diameter

## Angle measurement

mm

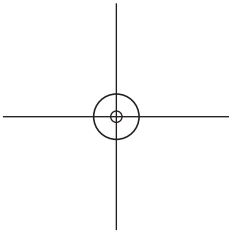
unit of measurement

## Additional Information

The circle is intended to explain the adjustment of magnification. The information given here refer to a PC monitor, not to the TFT display.

- For correct function of the software 2 things are to be considered:
  1. You must have administrator privileges on your computer, otherwise you get error messages.
  2. You must have installed PC DirectX9c on your computer. To determine whether DirectX9c is installed, enter the Start> Run the command „dxdiag“.
- The DigiMicro has two fixed focus settings. Turn the knurled screw accordingly. Near the right keystroke you get the low magnification, in the vicinity of the left keystroke you get high magnification.

Turn the thumbscrew to the right keystroke. Set a software resolution to 5M.

Set the DigiMicro Mobile with the illuminated front to the centre of the drawing (left).

Wait 2 seconds until the white balance is adjusted.

Turn the thumbscrew slowly to the left.

First you see the outer circle on full screen. This is the low magnification. It has a diameter of 6mm.

Now, if you continue turning the thumb wheel to the left, you see the inner circle full screen. This is the high magnification. It has a diameter of 1.5 mm.

### Magnifications

The information to the magnification refers exclusively to the representation on a monitor and not on prints.

The scale on the knurled screw serves for the input with the provided measuring software.

In praxis this ensures in a deviation of 2-5% max.

### Magnification 20x, 200x und 500x

20x:

The minimum magnification is near the right stop of the thumbscrew(rear view direction). In direct contact of the illuminated point of the microscope you will get a magnification of about 80-90x if the image is shown in fullscreen on a 17" monitor.

200x:

The minimum magnification is near the left stop of the thumbscrew(rear view direction). In direct contact of the illuminated point of the microscope you will get a magnification of about 290-300x if the image is shown in fullscreen on a 17" monitor.

500x:

The minimum magnification is near the left stop of the thumbscrew(rear view direction) and the

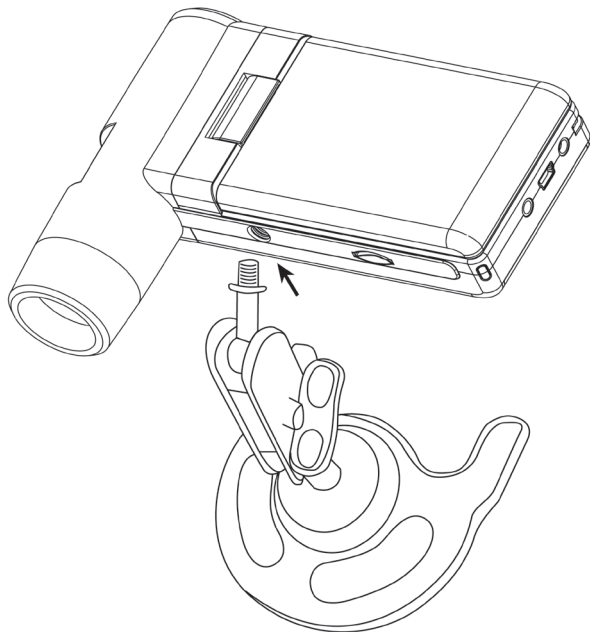
digital zoom of the DigiMicro Mobile is set to 2x-4x. In direct contact of the illuminated point of the microscope you will get a magnification of about 500-700x if the image is shown in fullscreen on a 17" monitor.

To get a fullscreen 100%(1:1) on the monitor please use a freeware program like XnView (xnView.de) or Irfanview (irfanview.de).

### **Mounting the stand**

Put the stand end with screw toward the mount hole on the microscope properly and rotate the end into the hold carefully. Do not force the end into the hole; forcing may damage the microscope.

Adjust the height of the stand at a proper position.

### **Service information**

In case of complaints your device shows a defect, please consult your dealer or contact the service department of Drahtlose Nachrichtentechnik Entwicklungs- und Vertriebs GmbH to agree a servicing.

Service time: Monday until Thursday from 8.00-17.00h  
Friday from 8.00-16.00h

Hotline/service: +49 6074 3714 31

E-mail: support@dtectr.de

Please do not send your device to our address without request by our service team. The expense and the risk of the loss are for debits of the sender. We reserve to refuse the unrequested shipment or return corresponding goods to the sender at buyer's risk and expense.

### Specifications

Image sensor	5 megapixels
Display	7,6 mm (3") TFT
Video resolution	12M(4032x3024), 9M (3456x2592), 5M(2592x1944), 3M(2048x1536), 1.3M(1280x960), VGA (640*480) Pixel
Still capture resolution	12M(4032x3024), 9M (3456x2592), 5M(2592x1944), 3M(2048x1536), 1.3M(1280x960), VGA (640*480) Pixel
Colour	24-Bit-RGB
Optic	Dual-Linse 27x + 100x
Focus range	manuel, 10 mm till 300 mm
Flicker Frequency	50/60 Hz
Frame rate:	max. 30 fps
Magnifications	ca 20 and 200x, 500x
Image format	JPG
Video format	AVI
White balance	automatic
Light source	8 LED, white with dimmer
PC interface	USB 2.0
Operating system	Windows XP SP2/Vista, Windows 7, Windows 8/10, MAC OS 10.6 or higher
Power supply	Li-Ionen battery 3,7V/800mA, 5 V (USB Port), charger
Dimensions	105 x 130 x 28 mm

### Declaration of Conformity

Hereby, Drahtlose Nachrichtentechnik Entwicklungs- und Vertriebs GmbH declares that this Digi-Micro Scale is in compliance with the essential requirements and other relevant provisions of Directive 2014/30/EU.

The declaration of conformity is available for download [www.dtectr.de/konformitaet](http://www.dtectr.de/konformitaet)

### Disposal

The article, its packaging and the batteries supplied have been manufactured from valuable materials that can be recycled. Recycling reduces the amount of waste and helps to protect the environment. Dispose of the packaging at a recycling point which sorts materials by type. Make use of the local facilities provided for collecting paper; cardboard and lightweight packaging.

You are legally obliged to dispose of old appliances separately from household refuse! Information about where they can be deposited free of charge is available from your local authorities.

Regular batteries and rechargeable batteries must not be disposed of along with household refuse! You are legally obliged to dispose of flat batteries or rechargeable batteries at a local authority collection point or to return them to a battery retailer.

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## **Warranty regulations**

1. In general the European warranty regulations are valid. Warranty starts with receipt of unit. In a warranty case, please submit the unit together with invoice, unit pass and error description. Without these documentary evidence we are not obliged to a warranty performance.
2. During warranty period we repair all constructional or material defects which are caused by a manufacturing fault. It is in the decision of the manufacturer to change or replace defective parts/units. Further claims are excluded. Changed parts will automatically go over into our property. We will not be held responsible for future damages.
3. Manipulation on the unit through non-authorized companies/persons automatically leads to an end of warranty period.
4. Excluded of warranty claims are damages due to improper treatment, not following the manual, oxidation of batteries, permanent usage.
5. The initial warranty period is not extended because of repair or replacement of parts.
6. These warranty regulations are valid in the Federal Republic of Germany.

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